

Getting started with Consibio Cloud

Consibio Cloud is your window into your Consibio devices: live measurements, historical data, alarms, reports and remote control — all in one place.

Sign in

1. Go to v3.consibio.cloud (<https://v3.consibio.cloud>).
2. Sign in with your e-mail address and password.
 - If your organization already uses Consibio Cloud, ask a project administrator to [invite you](#) — you'll receive an e-mail with a link to create your account.

Projects

Everything in Consibio Cloud lives inside a **project**: devices, sensor data, alarms, reports and team members. When you sign in you'll land in your default project. If you're a member of several projects, use the project switcher in the menu to change between them.

Find your way around

The main menu gives you access to:

- **Dashboard** — widgets with live and historical measurements from your devices.
- **Devices** — every device in the project, its status and battery level. See [Claiming a device](#).
- **Alarms** — rules that notify you when a measurement goes out of range. See [Creating an alarm](#).
- **Reports** — periodic summaries of your data.
- **Users** — the members of your project and their roles.
- **Settings** — project configuration.

Next steps

- [Claim your first device](#)
- [Create an alarm](#)
- [Invite your team](#)