

# Exporting data to CSV

Consibio Cloud can export your project's measurements as a CSV file, which opens directly in Excel or any other spreadsheet tool.

## Export data

1. Open the **Reports** view from the main menu.
2. Select **Export data**.
3. Under **Type**, keep **CSV export** selected.

### **Note**

PDF and Excel export types are planned but not available yet — CSV is currently the only option.

4. Under **Elements**, choose which measurements to include. If you don't select any, all elements are included.
5. Choose the **time range** the export should cover.
6. Optionally set **Align timestamps** to an interval (for example 60s) so every element's readings line up on the same rows. Leave it at 0s to export each element's raw, unaligned timestamps.
7. Select **Download**.

Your browser downloads a file named `Consibio Cloud Datalog.csv`.

## Opening the file in Excel

Double-click the downloaded file to open it in Excel. Depending on your Windows system language, Excel may not split the columns correctly on the first open — if that happens,

select column A, then use **Data > Text to Columns** in Excel's toolbar to split it, choosing **Comma** as the delimiter.

## Support

Need help with a data export or import? Contact [support@consibio.com](mailto:support@consibio.com) — we can help with imports on Mac, Windows, Linux and specific Excel versions.